Petite Digital Stirrer

English
Instruction Manual
pages 2-7

Espanõl
Manual de intrucciones
pages 14-19

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Contents

GENERAL INSTRUCTIONS.....	3
SAFETY INSTRUCTIONS	3
APPROVALS	4
TECHNICAL DATA.....	4
STIRRER DESCRIPTION.....	5
INSTALLATION.....	5
ENVIRONMENTAL OPERATING CONDITIONS.....	6
GENERAL OPERATING INSTRUCTIONS.....	6
CLEANING, SERVICE, TRANSPORT & STORAGE.....	6
WARRANTY AND PRODUCT LIABILITY.....	7

PACKING LIST

- Petite digital stirrer
- Power cord
- Stand base
- Stand support rod and lock nut
- Impeller shaft
- (2) Impeller blades
- (2) Impeller lock screws
- Instruction manual

250PK004
Rev0 - 01/08

GENERAL INSTRUCTIONS

Read all safety instructions before operating.

Unpack the stirrer carefully. Ensure the stirrer is not damaged and verify the contents against the packing list. If the stirrer is damaged or the contents do not correspond with the packing list, notify your supplier immediately.

Read the manual carefully! Make sure that all users read the manual thoroughly before operating the stirrer. For further questions or inquiries, please contact:

Caframo Ltd.

RR #2, Warton, Ontario, Canada N0H 2T0

Tel: (519) 534-1080

Toll free in USA and Canada: (800) 567-3556

Fax: (800) 209-6786

Email: labsales@caframo.com

Web site: www.caframo.com

SAFETY INSTRUCTIONS

Each operator should read all instruction before using this device.

The stirrer must be securely fixed to a stable support such as the provided stand. If other stands are used, the unit must be secured such that it will not fall if tipped.

CAUTION: This is not an explosion-proof stirrer. Do not use in a highly flammable or explosive environment.

Spinning paddles or impellers can cause severe personal injuries. Operators must use extreme care and good judgment when mixing at any speed.

All mixing paddles and impellers must be in good condition with straight shafts. If the stirrer vibrates at high speeds, check the paddle shaft for damage, and repair or replace if needed.

Extreme care must be taken when mixing chemicals to ensure that no chemicals are splashed outside the mixing vessel. Care must be taken when changing to faster mixing speeds.

Ensure that the mixing impeller does not contact the containment vessel.

Do not operate while standing in water. Keep the unit dry and do not immerse any part, except the mixing paddle into any liquids. Protect unit and operator from splashing.

Ensure that no loose clothing, jewelry, or hair can become entangled in any rotating parts. A fast spinning chuck can cause injury to operator.

Power can be interrupted to the stirrer by pressing the power button. In an emergency, disconnect the main power cord.

APPROVALS

North American

Constructed in accordance with CAN/CSA-C22.2 (1010),
FCC 15/ICES-003/EN 55011

Tested to the same standards as IEC61010.

European

Constructed in accordance with EMC/EMI standard EN61326-1:2006 (EN55011, conducted and radiated emissions, EN61000-3-2, EN61000-4-2, EN61000-4-3, EN61000-4-4, EN61000-4-5, EN61000-4-6)

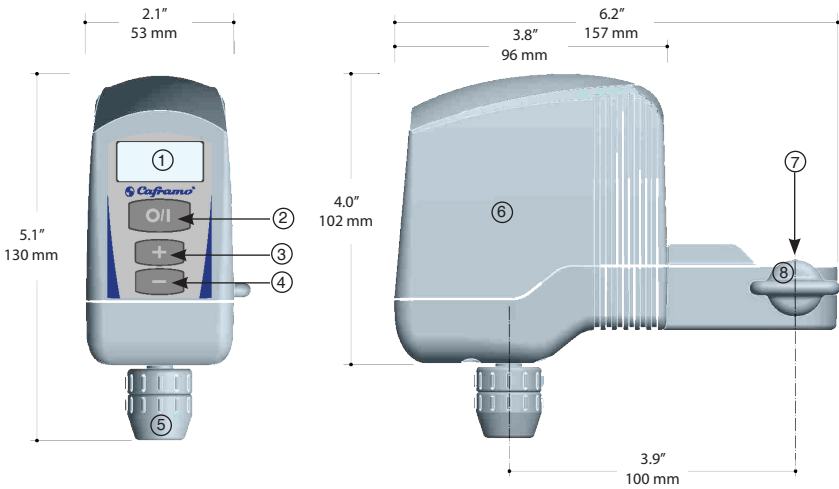
RoHS and WEEE compliant.

TECHNICAL DATA

Model Number	BDC250
Speed range	50 To 2500 rpm
Maximum torque	10 Ncm
Mechanical output power	1/100 hp, 7 watts
Input to power supply	AC 100 to 240 volts, 50/60 hz (must use supplied Caframo power cord)
Electrical output from power supply to stirrer	DC 24 V, 0.5 A
Weight - stirrer only	0.9 lbs (.4 kgs)
Weight - total package	3.1 lbs (1.42 kgs)
Maximum volume	2 liters, water-like viscosity
Molded housing	Chemical-resistant, high-impact nylon. Recyclable (PA).
Keypad	Waterproof, chemical-resistant polyester
Power cord length	6 ft
Keyless chuck	Fits 6 mm shaft
LED display accuracy	±10 rpm, 4 digit

Technical specifications subject to change without notice.

STIRRER DESCRIPTION

- | | | | |
|---|------------------------|---|---------------|
| ① | Four-digit LED display | ⑤ | Keyless chuck |
| ② | Button - power on/off | ⑥ | Housing |
| ③ | Button - increase rpm | ⑦ | Clamp opening |
| ④ | Button - decrease rpm | ⑧ | Clamp knob |

INSTALLATION

Stand Assembly

Remove lock nut. Insert the threaded end of the stand rod into the stand base. Turn rod clockwise to tighten. Secure with lock nut.

Stirrer Installation

Slide the clamp opening over the top of the stand rod. Lower to desired height. Turn clamp knob to securely tighten the stirrer onto the stand rod. Connect supplied power cord to stirrer. Connect power cord to a grounded electrical outlet.

Impeller Assembly

Add impeller blade(s) to impeller shaft. Tighten screw to secure.

Impeller Installation

Turn lower half of chuck to open. Do not remove completely. Insert impeller end into chuck opening until it stops. Turn lower half of chuck (finger tight) to close chuck and secure impeller.

ENVIRONMENTAL OPERATING CONDITIONS

The BDC250 stirrer must operate in the following conditions:

- Indoors
- Altitudes not exceeding 2000 meters
- Temperature from 41 to 104°F (5 to 40°C)
- Maximum relative humidity 80% for temperatures up to 87°F (31°C) decreasing linearly to 50% relative humidity at 104°F (40°C)
- Pollution degree 2 in accordance with IEC 664 “Normally only non-conductive pollution occurs. Occasionally, however, temporary conductivity caused by condensation may be expected.”

GENERAL OPERATING INSTRUCTIONS

The stirrer must be securely mounted as specified in the installation instructions. Ensure the mixing shaft and blade are securely attached. The stirrer is now ready for use to mix liquids.

Button Operation Functions:

Power Button - Turns the stirrer motor on and off.
When initially turned on, stirrer will spin at 50 rpm.

Increase Button - Increases the speed.

Decrease Button - Decreases the speed.

- OL Indicates an overload condition. The load is too great for the stirrer and must be reduced. To reset stirrer turn power off and back on.

CLEANING, SERVICE, TRANSPORT & STORAGE

Cleaning

Using a mild soap and water solution, wipe the housing and control panel of the stirrer with a damp cloth. Do not use chlorine bleach, chlorine-based cleanser, abrasives, ammonia, steel wool, or scouring pads with metal content or similar harsh solvents or abrasives. These may damage the surface of the instrument.

Service

Do not attempt to service or repair a Caframo stirrer. **If the stirrer housing is opened, the warranty becomes void.** Contact Caframo for return authorization and return instructions.

Transport and Storage

Clean stirrer so that it is free from any materials which may be harmful. Provide an MSDS or other safety data sheets for all chemicals used in conjunction with this stirrer where appropriate. Place the stirrer and its parts into the original packing to protect and prevent damage during transport. Seal the original packing with packing tape. Store the packed stirrer in a dry place.

CAUTION: Failure to clean, service, and handle the stirrer as outlined above can cause it to fail which could impair the safety of the user.

WARRANTY AND PRODUCT LIABILITY

Warranty

Caframo warrants the herein described product of its manufacture for three years from receipt of the equipment by the purchaser, against defects in material and workmanship. Warranty registration is available on-line at www.caframo.com or fill out the supplied warranty registration card.

This limited warranty covers parts and labor. **If the stirrer housing is opened, the warranty becomes void.**

This warranty may only be altered by a specifically published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Expendable items are not covered by this warranty. Damages due to corrosion or accidental breakage are not covered.

Liability

Under no circumstances shall Caframo be liable for indirect, consequential, or special damages of any kind.

Caframo reserves the right to make technical changes without prior notice.

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